CHECKIT® Comparator

Chlor frei, gebunden, gesamt 0 - 3.5 mg/l

Chlorine free, combined, total 0 - 3.5 mg/l

Chlore libre, combiné, total 0 - 3.5 mg/l

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



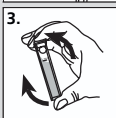
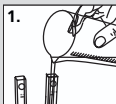
Note:

Read the safety data sheets for the chemicals used. Use the recommended personal protective equipment and follow the disposal instructions.

Place the CHECKIT® Disc in the CHECKIT® Comparator with numeric values facing the operator.

**VARIO Free Chlorine**

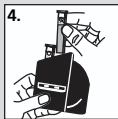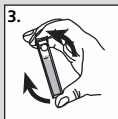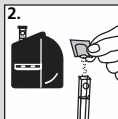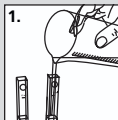1. Rinse the cuvettes with the water sample. Then fill both cuvettes with the water sample up to the 5 mL mark. Close one cuvette with the lid and place it in the left chamber of the CHECKIT® Comparator as a blank sample. The dot on the cuvette must point to the viewer.
2. Add the powder from the VARIO DPD Free Chlorine F5 packet to the other cuvette.
3. Close the cuvette with the cap and gently swirl to dissolve the powder.
4. Place the cuvette in the right chamber of the CHECKIT® Comparator. The dot on the cuvette must point to the viewer.
5. Match the colour of the water sample with the CHECKIT® DISC. The colour corresponding to the water sample indicates the free chlorine content in mg/L.

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1. Rinse the cuvettes with the water sample. Then fill both cuvettes with the water sample up to the 5 mL mark. Close one cuvette with the lid and place it in the left chamber of the CHECKIT® Comparator as a blank sample. The dot on the cuvette must point to the viewer.
2. Add the powder from the VARIO DPD Total Chlorine F5 packet to the other cuvette.
3. Close the cuvette with the cap and gently swirl to dissolve the powder.
4. Wait for the 3 minutes reaction time.
5. Place the cuvette in the right chamber of the CHECKIT® Comparator. The dot on the cuvette must point to the viewer.
6. Match the colour of the water sample with the CHECKIT® DISC. The color corresponding to the water sample indicates the total chlorine content in mg/L.

Combined Chlorine

Total chlorine value (mg/L) – free chlorine value (mg/L) = combined chlorine value (mg/L).

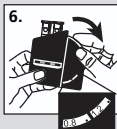
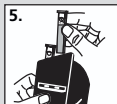
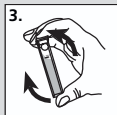
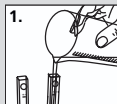
Note:

Levels of chlorine above 10 mg/L can cause bleaching of the sample.

Important:

After each measurement, rinse the cuvettes thoroughly. Store CHECKIT® Disc in the dark.

If stored properly, the color discs have a shelf life of 3 years from the date of receipt.

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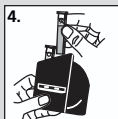
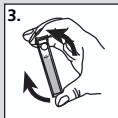
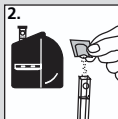
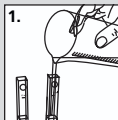
Chlore combiné

Concentration de chlore total (mg/L) - Concentration de chlore libre (mg/L) = Concentration de chlore combiné (mg/L)

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