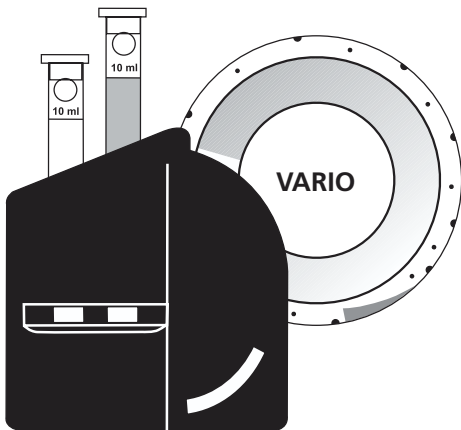
CHECKIT® Comparator

Mangan LR 0.1-0.7 mg/L Mn

Manganese LR 0.1-0.7 mg/L Mn

Manganèse LR 0.1-0.7 mg/L Mn

Manganeso LR 0.1-0.7 mg/L Mn

DE

EN

FR

ES

IT

Hinweis:

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




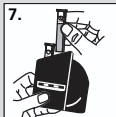
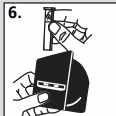
Note:

Read the safety data sheets for the chemicals used. Use the recommended personal protective equipment and follow the disposal instructions.

Place the CHECKIT® Disc in the CHECKIT® Comparator with numeric values facing the operator.

**VARIO Manganese**

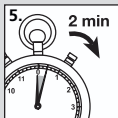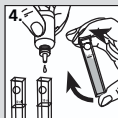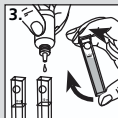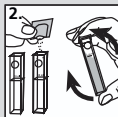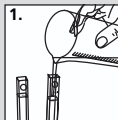1. Fill one cuvette with the water sample up to the 10 mL mark. Fill the other cuvette with DI water up to the 10 mL mark.
2. Add one powder from the VARIO Asorbic Acid packet to each of the two cuvettes. Close the cuvettes with the caps and gently swirl to dissolve the powder.
3. Add 15 drops Alkaline-Cyanide Reagent Solution to each of the two cuvettes. Hold the bottle vertically. Close the cuvettes with the caps and gently swirl to mix.
4. Add 21 drops PAN Indicator Solution to each of the two cuvettes. Hold the bottle vertically. Close the cuvettes with the caps and gently swirl to mix.
5. Wait for the 2 minutes reaction time.
6. Place the cuvette with DI water in the left chamber of the CHECKIT® Comparator as a blank sample. The dot on the cuvette must point to the viewer.



7. Place the cuvette with water sample in the right chamber of the CHECKIT® Comparator. The dot on the cuvette must point to the viewer.
8. Match the colour of the water sample with the CHECKIT® Disc. The colour that matches the water sample indicates the Manganese content in mg/L.

Note:

For water samples with a hardness of more than 300 mg/L, add 10 drops of VARIO Rochelle Salt Solution after adding the powder. Then continue with step 3.

Important:

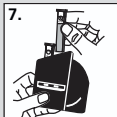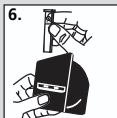After each measurement, rinse the cuvettes thoroughly. Store CHECKIT® Disc in the dark. If stored properly, the color discs have a shelf life of 3 years from the date of receipt.

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